



QuikRead[®] Connectivity

Improve traceability and eliminate data recording errors

- Connectivity to Laboratory/Hospital Information System
- Patient identification with Barcode reader
- Printing of patient results

POC

Take full advantage of your QuikRead® 101!

The QuikRead point-of-care system consists of a small photometer and assay-specific reagent kits. The test results are available within minutes. QuikRead Connectivity increases patient security and result traceability by eliminating risks involved in manual recording of test results.

Product	Cat. No.
QuikRead 101 Instrument	06040
QuikRead QuikLink Setup Wizard	06100
QuikRead QuikLink bar code reader (Datalogic DLC7070-M1)	06109



QuikRead 101 + pc

Transfer the QuikRead test result to your pc using a terminal program (for example HyperTerminal). This way you can save the test result in a text file and print daily summaries of the accrued data. Using a barcode reader and QuikRead QuikLink, test results can be coupled with patient identifications.




QuikRead 101 + printer

Connect your QuikRead to a printer to obtain a printout of the test result. Manual data recording errors can be eliminated by attaching the printout to the patient's medical record.



QuikRead 101 + barcode reader

Using a barcode reader and QuikRead QuikLink, test results can be coupled with patient identifications.

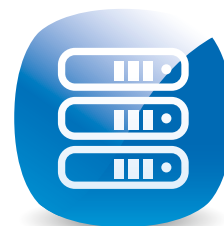

QuikRead 101 + LIS/HIS

Integrate your QuikRead with an electronic patient record or a laboratory/hospital information system (LIS/HIS) using QuikRead QuikLink. QuikRead QuikLink allows test results coupled with patient identifications to be transferred from a point-of-care site to a central database.



QuikRead 101 QuikLink

QuikRead QuikLink consists of QuikRead QuikLink Setup Wizard and a separate barcode reader. Your QuikRead 101 Instrument can be configured with QuikRead QuikLink to communicate in a standard format with an electronic patient record or LIS/HIS.

